

URVALUE	GENERAL LEVEL	Lab Administrator	Lab Analyst	Lab Supervisor	Administrative Assistant	Lab Administrator	EXPANDED DESCRIPTION
ACCPSTRT APP	ADMIN / USER	X					Allows the user to run the Access Report program
ADD_REPLICATE	ADMIN / USER	X	X	X	X		Allow the user to add new Replicate
AER_RUN	ADMIN	X					Allows the system to run automated email reports
AMAIN1 APP	ADMIN / USER	X	X	X	X		Allows the user to run the Analysis List Maintenance Program from the LABWORKS Desktop
ANALVALIDATE	ADMIN / USER	X		X			Allows the user to validate analysis
ANALYSIS_STATE	ADMIN	X					Allows the user to modify the analysis state in Analysis List Maintenance
ANLOBJECTCHANGE	ADMIN / USER	X	X	X			Allows the user to modify or delete an object stored with an analysis
ANLOBJECTLOAD	ADMIN / USER	X		X			Allows the user to load an object onto an analysis
ANLOBJECTVIEW	ADMIN / USER	X	X	X			Allows the user to view an object that has been stored on an analysis
APPROVE PRIO	ADMIN / USER	X		X	X		Run Object Publisher
ARCHIVE APP	ADMIN	X					Allows the user to run the Sample Archive Program from the LABWORKS Desktop
ASSIGN ANALYST	ADMIN / USER	X	X	X			Allows the user to assign an analyst to an analysis
ASSIGN DUEDATE	ADMIN / USER	X		X			Allows the user to assign a due-date on an analysis
ASSIGN INSTR	ADMIN / USER	X		X			Allows the user to assign an instrument to an analysis
AUDIT_CHKSUMVIEW	ADMIN	X			X		Allows the user to view the Check-sum field within the Audit-trail viewer
AUDIT_CHKSUM	ADMIN	X					Allows the user to update the check-sum field on an audit record
AUDIT_PIV	ADMIN / USER	X					Allows the user to access the Audit-trail Viewer
AUDITRUN	ADMIN / USER	X	X	X	X		Allows the user to run the Audit-trail viewer
BACKLOG APP	ADMIN / USER	X	X	X			Allows the user to run the Backlog Program from the LABWORKS Desktop
BACKLOGRUN	ADMIN / USER	X		X			Allows the user to run a backlog report
BARCODERUN	ADMIN / USER	X	X	X			Allows the user to generate a barcode (within the Sample/Modify Program)
BATCHDELETE	ADMIN	X					Delete QA/QC Batches
BATCHSETUP	ADMIN / USER	X	X				Create QA/QC Batches
CALCSETUP	ADMIN	X					Allows the user to setup / create calculations in the lins
CAL_MODIFY	ADMIN / USER	X					Allows the user to modify CAPA Reports
CHANGEPASSWORD	ADMIN / USER	X	X	X	X		Allows the user to change their password
CINV	ADMIN / USER	X					Allows the user to run the Chemical Inventory program
CINV_APPR ITEMS	ADMIN / USER	X					Allows user to Approve Item Stock in Chemical Inventory
CINV_CR_ITEMNAME	ADMIN / USER	X					Allows user to create new item in Chemical Inventory
CINV_CR_VENDOR	ADMIN / USER	X					Allows user to Create Vendor
CINV_MODIFY ITEM	ADMIN / USER	X					Allows user to edit Chemical Inventory Item Stock
CINV_OVRDEXPDATE	ADMIN / USER	X					Allows user to override expiry date of item in Chemical Inventory
CINV_PR_APPR	ADMIN / USER	X					Allows user to Approve Purchase Requisition
CINV_RECONCILE	ADMIN / USER	X					Allows user to Physical Reconciliation
CINV_SPUSER	ADMIN / USER	X					Chemical Inventory Super User
COA BYPASS VIOLS	ADMIN / USER	X					Allows the user to generate a COA that has specification violations
COADLETE	ADMIN	X					Allows the user to delete a COA Customer Base Record
COAPROD APP	ADMIN / USER	X					Allows the user to run the COA Program from the LABWORKS Desktop
COARUN	ADMIN / USER	X					Allows the user to run / generate a COA
COASETUP	ADMIN	X					Allows the user to configure the COA program
CONFIGDESKMENU	ADMIN	X		X			Allows the user to configure the LABWORKS Desktop Menu
CONFIGWEBMENU	ADMIN	X		X			Allows the user to configure the LABWORKS Webtool Menu
CONVERTCALCDS	ADMIN	X					Allows the user to run the ConvertCalc Program (update calculations from 6.3 to 6.4)
CONVSPEC00T01	ADMIN	X					Allows the user to run the ConvertSpecs Program (update specifications from 6.3 to 6.4)
CRPSET APP	ADMIN	X					Allows the user to run the Crystal Report Setup program from the LABWORKS Desktop
CTRACK APP	ADMIN / USER	X	X	X	X		Allows the user to run the Custody Tracking program from the LABWORKS Desktop
CUSTODYRUN	ADMIN / USER	X	X	X	X		Allows the user to enter Custody Tracking information into Custody Tracking
DELETE_REPLICATE	ADMIN / USER	X	X	X	X		Allow the user to delete existing Replicate
	ADMIN	X					Allows the user to register and modify the registration of programs associated with the Labworks Desktop (pre 6 privilege)
DESKTOPSETUP	ADMIN	X					
DMR APP	ADMIN / USER	X					Allows the user to run the DMR application from within the LABWORKS Desktop
DMRREAD	ADMIN / USER	X					Allows the user to view DMR information in the DMR Application
DMRRUN	ADMIN / USER	X					Allows the user to create DMR reports information in the DMR Application
DMRSETUP	ADMIN	X					Allows the user to setup DMR information in the DMR Application
DMRWRITE	ADMIN / USER	X					Allows the user to enter DMR information in the DMR Application
DOCKCREATE	ADMIN	X					Allows the user to create new documents in the Documentation tracking program
DOCKDELETE	ADMIN / USER	X					Allows user to delete Document Reference
DOCKINFO APP	ADMIN / USER	X	X	X	X		Allows the user to run the Documentation tracking program from the LABWORKS Desktop
DOCKREFVIEW	ADMIN / USER	X					Allows the user to review Documents
DOCKVIEW	ADMIN / USER	X	X	X	X		Allows the user to view documents referenced in the Documentation tracking program
DV_SHOW_HISTORY	ADMIN / USER	X	X	X	X		Allows the user to Show History from the Dataviewer
ERRORLOGDELETE	ADMIN	X					Allows the user to delete errorlog entries
ERSETUP APP	ADMIN / USER	X	X	X	X		Allows the user to run the Execution Report setup program from the LABWORKS Desktop
EWKSHKTCOMPLET	ADMIN / USER	X					Allows the user to complete eWorksheet
EWKSHKDELETE	ADMIN / USER	X					Allows user to delete eWorksheet
EWKSHKEDIT	ADMIN / USER	X					Allows the user to Change eWorksheet to In Progress
EWKSHKSETUP	ADMIN / USER	X					Allows user to setup eWorksheet
EWKSHKTEMPXCEL	ADMIN / USER	X					Allows user to Edit eWorksheet excel template
EWKSHKTEMPSETUP	ADMIN / USER	X					Allows user to setup eWorksheet template
EWKSHKVALIDATE	ADMIN / USER	X					Allows the user to Validate eWorksheet
EXCEPTIONRUN	ADMIN / USER	X	X	X	X		Allows the user to run Exception Reports from the Exception Report program
	ADMIN	X					Allows the user to configure the Automatic Exception Reporting program. This program will detect exceptions and report them with a stored Word document, email or sms the exceptions to a user or a group of users
EXCPTRIPROTESTSETUP	ADMIN / USER	X					Allows the user to run the Automatic Exception Report program from the LABWORKS Desktop
EXCPTRLO APP	ADMIN / USER	X	X	X	X		Allows the user to run the Export Data program from the LABWORKS Desktop
EXPORT APP	ADMIN / USER	X	X	X	X		Allows the user to export data from the LIMS using previously configured export formats
EXPORTRUN	ADMIN / USER	X	X	X	X		Allows the user to export data from the LIMS into some required format. Several options for exporting are available. See the User Manual.
EXPORTSETUP	ADMIN	X					Allows the user to run the Field Entry Data Terminal program from the LABWORKS Desktop
FDSET APP	ADMIN / USER	X					Allows the user to print labels to be used for field sampling.
FIELDLABELRUN	ADMIN	X					Allows the user to create special lab forms using the Forms Generator Program
FORMSETUP	ADMIN	X				X	Allows the user to configure Persisted Group information in LWSampleLogin
GROUPTYPECONFIG	ADMIN	X					
IMANAGE APP	ADMIN / USER	X	X	X			Allows the user to run the Instrument Maintenance and Calibration program from the LABWORKS Desktop
IMCHODIV	ADMIN	X		X			Allows the user to modify the data within the Instrument and calibration database
INEDIT	ADMIN	X			X		Allows the user to change the INI (initialization) file
INSTCONV APP	ADMIN	X					Allows the user to run the Instrument conversion program from the LABWORKS Desktop
INSTUB APP	ADMIN	X					Allows the user to run the Instrument Library program from the LABWORKS Desktop
INVOICERUN	ADMIN / USER	X	X	X	X		Allows the user to run / generate invoices using the standard Labworks Invoice program.
	ADMIN / USER	X					Allows the user to run the Industrial Pretreatment Program. (IPP) This is a US EPA regulatory program that tracks the permits of industrial pretreatment into a Publicly Owned Wastewater Treatment Facility
IPPRUN	ADMIN / USER	X					
IPPSSETUP	ADMIN	X					Allows the user to configure the IPP program
LIPO APP	ADMIN / USER	X	X	X			Allows the user to run the Labworks Industrial Pretreatment Option
LIANINT APP	ADMIN / USER	X	X	X			Allows the user to run the Location List Maintenance program from the LABWORKS Desktop
LOCATIONMODIFY	ADMIN	X	X	X			Allows the user to setup the location codes in the Location List Maintenance program.
LOCATIONHREAD	ADMIN / USER	X	X	X			Allows the user to view / print location code information from Location List Maintenance
LOGINWOLDCODE	ADMIN	X					Allows the user to login samples that do not have a predefined Location Code
LOGPRINTSETUP	ADMIN	X					Allows the user to configure the information that may be printed a sample login
LOGTEMPPLATEDEL	ADMIN	X					Allows the user to delete a Sample Login Template
LOGTEMPPLATESETUP	ADMIN / USER	X	X	X	X		Allows the user to configure the single sample, multi sample and Process/Scheduler sample login screens or templates
LWEXP_INTERNAL	ADMIN / USER	X	X	X	X		Labworks Explorer - Users without this priv don't see samples in LWExplorer until they are Validated.
LWEXPCONFIGURE	ADMIN / USER	X	X	X	X		Allows the user to configure LABWORKS Explorer
LWEXPRUN	ADMIN / USER	X	X	X	X		Allows the user to run LABWORKS Explorer
LWEXPSHOWRAW	ADMIN / USER	X	X	X			Allows the user to view un-validated data using the read only Labworks Explorer program
LWSEQC APP	ADMIN / USER	X	X	X	X		Allows the user to run the LWSEQC program from the LABWORKS Desktop
LWTFREND APP	ADMIN / USER	X	X	X	X		Allows the user to run the LWTFREND program from the LABWORKS Desktop
LWUPDATES	ADMIN	X			X		Allows the user to apply database updates in the LWMeritonUtility program
LWWEDEF APP	ADMIN / USER	X					Allows the user to run the DEF program from the LABWORKS

LWWRMDE APP	ADMIN / USER	X					Allows the user to run the Form Editor program from the LABWORKS Desktop
LWWINDE APP	ADMIN	X				X	Allows the user to run the INI Editor program from the LABWORKS Desktop
LWWINV APP	ADMIN / USER	X	X	X	X		Allows the user to run the Invoice program from the LABWORKS Desktop
LWWMULTI APP	ADMIN / USER	X	X	X	X		Allows the user to run the Multi-sample Login program from the LABWORKS Desktop
LWWSMGL APP	ADMIN / USER	X	X	X	X		Allows the user to run the Single-sample login program from the LABWORKS Desktop
LWWSYSTA APP	ADMIN	X	X	X	X		Allows the user to run the System Manager program from the LABWORKS Desktop
MAILLISTSETUP	ADMIN	X	X	X	X		Allows the user to setup mail list records in SampleMailist
MAKETEMPLAT	ADMIN	X				X	Allows the user to create new templates under Demographic Data Entry
MCSERVO APP	ADMIN / USER	X	X	X			Allows the user to run the multi-component transfer program from the LABWORKS Desktop
METHODREFIDT	ADMIN / USER	X					Allows user to add custom Method Reference.
MIGRATE DATA	ADMIN / USER	X					Allows the user to Run migration tool application
MISSETUP	ADMIN	X					Allows the user to setup special files for calibration/askie mixture files
MODIFY REPLICATE	ADMIN / USER	X	X	X	X		Allow the user to change Replicate policy
MOR RPT APP	ADMIN / USER	X	X	X			Allows the user to run the MOR Report program from the LABWORKS Desktop
MORRUN	ADMIN / USER	X	X	X			Allows the user to run / generate MOR Reports from the MOR Report program
MORSETUP	ADMIN	X	X	X			Allows the user to configure the MOR (Excel based reporting tool)
MORSTCLE APP	ADMIN	X	X	X			Allows the user to run the MOR Setup program from the LABWORKS Desktop
MSAMPLES APP	ADMIN / USER	X	X	X			Allows the user to run the Modify/Delete Samples program from the LABWORKS Desktop
NGICREATEMODIFY	ADMIN	X					Allows the user to modify/create NGII Instrument Interfaces
NULL REPL VALUE	ADMIN / USER	X					Allows user to post Null Replicate result value
PAMRUN	ADMIN	X					Allows the user to use the Project Analytic Management (PAM) program to specify reortine limits on samples
PAMSETUP	ADMIN	X					Allows the user to configure the PAM program
PANAGATO APP	ADMIN / USER	X					Allows the user to run the Project Analytic Management Program
POETSETUP	ADMIN / USER	X					Allows the user to use the Portable Data Entry program (PDET)
PG PRIV	ADMIN	X					Allows the user to access the Persisted Groups in LWSampleLogin
PICKLISTSETUP	ADMIN	X					Allows the user to create exsists for use in certain Labworks programs and forms
PICKLIST APP	ADMIN / USER	X	X	X	X		Allows the user to run the Picklist Maintenance program from the LABWORKS Desktop
POST REPL VIOL	ADMIN / USER	X					Allows user to test Replicate results with violation
PRELIDC MODIFY	ADMIN	X					Allows the user to modify pre login group information
PREPORT APP	ADMIN / USER	X	X	X	X		Allows the user to run the Proares Report program from the LABWORKS Desktop
PRICULSTDELETE	ADMIN	X					Allows the user to delete a price list in the Location List Maintenance program
PRICULSTSETUP	ADMIN	X					Allows the user to create special prices lists for use in the Invoice program. Linked to special clients and customers so that discounts and or special billing can be managed.
PRINT FROM MOD	ADMIN / USER	X	X	X			Allows the user to print routing and work sheets from the modify and delete program
PROGRESSRUN	ADMIN / USER	X	X	X	X		Allows the user to run a Progress Report from the Progress Report program
PROGRESSSETUP	ADMIN	X	X	X			Allows the user to configure the Progress Report templates for customized viewing of the Labworks data
PROJECT MODIFY	ADMIN	X					Allows the user to modify Projects
PROJECT DELETE	ADMIN	X					Allows the user to delete Projects
PSEDTSCHEDULE	ADMIN	X	X	X			Allows the user to configure the Process Scheduler scheduling tool to create an automated sampleline plan.
PSLABRCFPT	ADMIN / USER	X	X	X	X		Allows the user to acknowledge Laboratory Receipt of samples in Process Scheduler
PSPRELADDELETE	ADMIN / USER	X	X	X	X		Allows the user to delete samples in Process Scheduler that do not have a LABWORKS Sample number assigned
PSPRELABMODIFY	ADMIN / USER	X	X	X	X		Allows the user to modify samples in Process Scheduler before a LABWORKS Sample number is assigned
PSPRODUCTOK	ADMIN / USER	X	X	X	X		Allows the user to acknowledge Production OK of samples in Process Scheduler
PSREORDER	ADMIN / USER	X					Allows the user to re-order samples within Process Scheduler
PSRESET COLUMN	ADMIN / USER	X					Allows the user to reset the column configuration within Process Scheduler
PSSAMPLELOGIN	ADMIN / USER	X	X	X	X		Allows the user to login samples within Process Scheduler
PSSHLELUFDEL	ADMIN	X					Allows the user to delete Shelf-life groups within Process Scheduler
PSSHLELUFEDIT	ADMIN	X					Allows the user to edit Shelf-life groups within Process Scheduler
PSSHLELUFEMAK	ADMIN / USER	X					Allows the user to make Shelf-life groups within Process Scheduler
PTMODIFY	ADMIN	X					Allows the user to modify the data within the Personnel Training database
QABATCH APP	ADMIN / USER	X	X	X			Allows the user to run the QA Batch Builder within the LABWORKS Desktop
QACRCRPO APP	ADMIN / USER	X	X	X			Allows the user to run the QACRC Report program from the LABWORKS Desktop
QATEMPL APP	ADMIN	X	X	X			Allows the user to run the QA Template Setup program from the LABWORKS Desktop
QATEMPLSETUP	ADMIN	X	X	X			Allows the user to setup QA Templates from the QA Template Setup program
QE DELETEQUERY	ADMIN	X					Allows the user to delete previously saved Query configured using the Query Engine
QE_EDITDATAMODEL	ADMIN	X					Allows the user to modify an existing Data Model (collection of tables containing the data being searched and the relationships required to allow that data to be retrieved accurately). This requires a person to understand the data structure and to have some knowledge of relational databases.
QE_EDITQUERY	ADMIN	X					Allows the user to modify an existing Query to change the output of the query
QE_NEWDATAMODEL	ADMIN	X					Allows the user to create a new data model by selecting the tables required to populate the queries that will be designed to extract data from the LIMS. This requires knowledge of the data structure and a basic understanding of relational databases.
QE_NEWQUERY	ADMIN	X					Allows the user to create a new query based upon an existing Data Model.
QE_RUNQUERY	ADMIN / USER	X					Allows the user to run an existing query from the Query Engine
QENGINE APP	ADMIN / USER	X					Allows the user to run the Query Engine program from the LABWORKS Desktop
RENTRY APP	ADMIN / USER	X	X	X	X		Allows the user to run the Result Entry program from the LABWORKS Desktop
REPORTRUN	ADMIN / USER	X	X	X	X		Allows the user to generate final reports using Labworks standard reporting tools
REPORTSETUP	ADMIN	X					LEGACY PRIVILEGE THAT DOES NOT USE CURRENT REPORTING TOOLS
RESET CONIG	ADMIN	X					Allows the user to reset configuration of all users
RESULTCHANGE	ADMIN / USER	X	X	X			Allows the user to modify a result value that has been saved into the database.
RESULTENTER	ADMIN / USER	X	X	X			Allows the user to load data, import data and calculate results. Required for almost all Labworks users who can add results to a sample
RESULTMERGE	ADMIN / USER	X					Allows user to merge previously posted result
RLTTEMPLATESETUP	ADMIN	X					Allows the user to create, modify and configure new or existing result entry templates to meet a users need for a custom data entry screen
RPT_TMPLT STATE	ADMIN / USER	X	X	X			Allows the user to update the Report Template State
RSETUP APP	ADMIN / USER	X	X	X			Allows the user to run the Result Entry Template program from the LABWORKS Desktop
SAMPARCHIVE	ADMIN	X					Allows the user to archive data using the Labworks data archival program
SAMPCOLLECTGRP	ADMIN / USER	X					Allows user to create or receive sample collection groups
SAMPDELETE	ADMIN	X					Allows the user to delete samples from the database. Once deleted, there is no trail for that sample. The audit trail ends with the sample below "deleted"
SAMPINACTIVATE	ADMIN	X	X	X	X		Allows the user to set a sample to an inactive status
SAMPLOCKOVERRIDE	ADMIN	X	X				Allows the user to modify stored sample and result information if the LOCKVALIDSAMPLES System Manager key is turned on.
SAMPLOGIN	ADMIN / USER	X	X	X	X		Allows the user to create samples in the LIMS. This is not the same as adding a sample at "not ready to collect" or at "ready to collect" status in Process Scheduler
SAMPMODIFY	ADMIN / USER	X	X	X			Allows the user to modify a sample after it has been loaded into the system
SAMPRESTORE	ADMIN	X	X	X			Allows the user to restore a sample to the active database using the Labworks Restore program
SAMPVALIDATE	ADMIN / USER	X	X				Allows the user to approve the data prior to making it available for reortine or for a Certificate of Analysis
SAVEDSKYTOP	ADMIN	X					LEGACY PRIVILEGE THAT IS NOT CURRENTLY USED
SAVEDSKYCREATE	ADMIN	X					CUSTOM Program for a customer, no general applicability
SAVEDSKYVIEW	ADMIN	X					CUSTOM Program for a customer, no general applicability
SCANDAEON APP	ADMIN	X	X	X			Allows the user to run the ScanDaemon program from the LABWORKS Desktop
SDGROUPSETUP	ADMIN	X					Allows the user to create and modify ScanDaemon Groups
SEND EMAIL	ADMIN	X				X	Allows the user to send an email through the LABWORKS Desktop
SEND MESSAGE	ADMIN / USER	X	X	X	X		Allows the user to send a message through the LABWORKS Desktop
SET EMAIL SERVER	ADMIN	X					Allows the user to configure email settings in the LABWORKS Desktop
SI SETUP APP	ADMIN	X					Allows the user to run the Special Information Form Designer from the LABWORKS Desktop
SIMLENTRYMODIFY	ADMIN	X					Allows the user to modify records in the Sample Mail List Entry
SPECIALINFOSETUP	ADMIN	X					Allows the user to configure a Special Information Form (SIFO) which can be used at the sample level and at the analysis code level
SPECSSETUP	ADMIN	X					Allows the user to set the specifications for location code analysis codes. Location code only to be sample specific or at the Primary Analysis code level to be a global specification for all samples using that analysis code
SQ CONFIGURE	ADMIN / USER	X	X	X			Allows the user to configure Stopped Queries
SO RUN	ADMIN / USER	X					Allows user to run Stopped queries
SO SAVE PRIVATE	ADMIN / USER	X	X	X			Allows the user to save 'Stopped Queries' to their Private profile
SO SAVE PUBLIC	ADMIN / USER	X	X	X			Allows the user to save 'Stopped Queries' for use by LABWORKS User
SQDRUN	ADMIN / USER	X	X	X	X		Allows the user to generate SQC Charts or use the SQC program as configured.
SQSETUP	ADMIN	X	X	X			Allows the user to configure the SQC Charting program for all normal Labworks SQC configuration attributes.
STATUSRUN	ADMIN	X					Allows the user to generate the system status report. A very general summary of the LIMS sample volume
STORRESULTSETUP	ADMIN	X					Allows the user to modify the system result abbreviations list
SYS CONFIGMENU	ADMIN	X				X	Allows the user to modify the System Menu's in the LABWORKS Desktop
TESTMODIFY	ADMIN	X	X	X			Allows the user to modify the definition of analysis codes that are previously configured in the database
TESTREAD	ADMIN / USER	X	X	X	X		Allows the user to view or print the details of an analysis code
TRAINING APP	ADMIN / USER	X					Allows the user to run the Personnel Training program from the LABWORKS Desktop