
LABWORKS LIMS v6.8

How to create Scheduled Task for CRGEN

Last Updated: March 30, 2018



How to create a Scheduled Task for CRGEN

1. Create and Test a BAT file to create batch file with CRGEN parameters.


Example:

```
"C:\LABWORKS\CLIENT\CRGEN6.exe" USER:USR PW:1 DB:DEFAULT  
INI:C:\LABWORKS\LWDATA\CRYSTAL\CR_SIMPLE.INI SL:c:\labworks\lwuser\x.dat
```

In this example the sample list is from a sample list previously created.

2. Confirm the batch file runs when the Labworks desktop isn't running, and the report is generated without prompts.

Create Basic Task

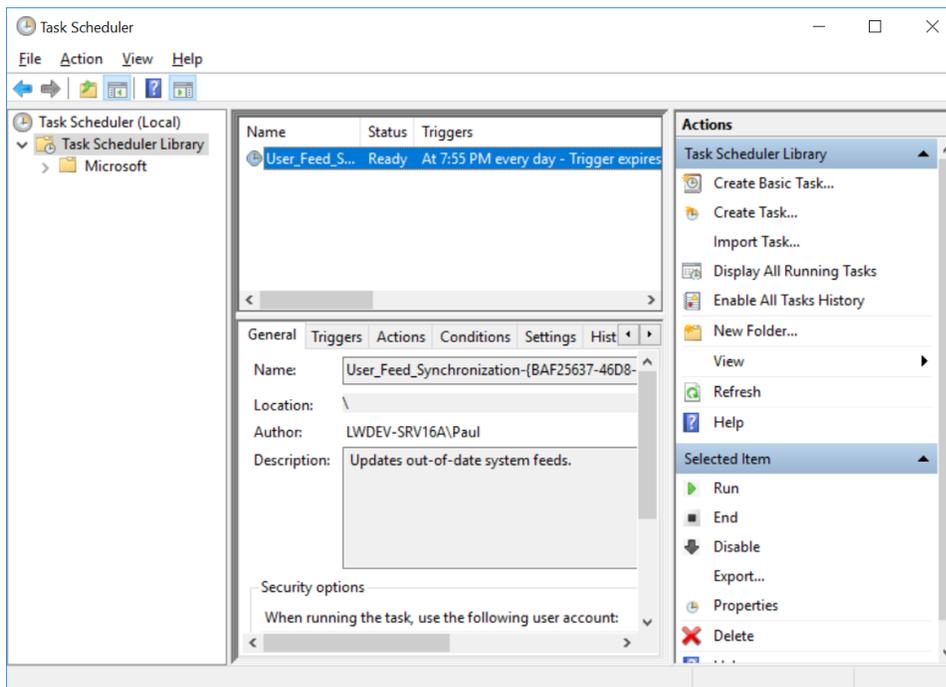

Enter Name/Description

Create Basic Task Wizard

Create a Basic Task

Use this wizard to quickly schedule a common task. For more advanced options or settings such as multiple task actions or triggers, use the Create Task command in the Actions pane.

Create a Basic Task

Trigger

Action

Finish

Name: CRGEN

Description: CRGEN Report

< Back Next > Cancel

Define the Schedule

Create Basic Task Wizard

Task Trigger

When do you want the task to start?

Trigger

Daily

Action

Finish

Daily

Weekly

Monthly

One time

When the computer starts

When I log on

When a specific event is logged

< Back Next > Cancel

Schedule Continued...

The screenshot shows the 'Create Basic Task Wizard' dialog box at the 'Daily' step. The window title is 'Create Basic Task Wizard'. On the left, there is a vertical list of steps: 'Trigger' (selected), 'Action', and 'Finish'. The 'Daily' step is highlighted with a blue bar. The main area contains the following fields and options:

- Start:** 11/17/2017 (calendar icon) 3:30:00 PM (time dropdown)
- Synchronize across time zones
- Recur every:** 1 days

At the bottom right, there are three buttons: '< Back', 'Next >', and 'Cancel'. The 'Next >' button is highlighted with a blue border.


Select Program to Run

The screenshot shows the 'Create Basic Task Wizard' dialog box at the 'Action' step. The window title is 'Create Basic Task Wizard'. On the left, there is a vertical list of steps: 'Trigger', 'Action' (selected), and 'Finish'. The 'Action' step is highlighted with a blue bar. The main area contains the following options:


- What action do you want the task to perform?**
- Start a program
- Send an e-mail (deprecated)
- Display a message (deprecated)

At the bottom right, there are three buttons: '< Back', 'Next >', and 'Cancel'. The 'Next >' button is highlighted with a blue border.


Select program to run...

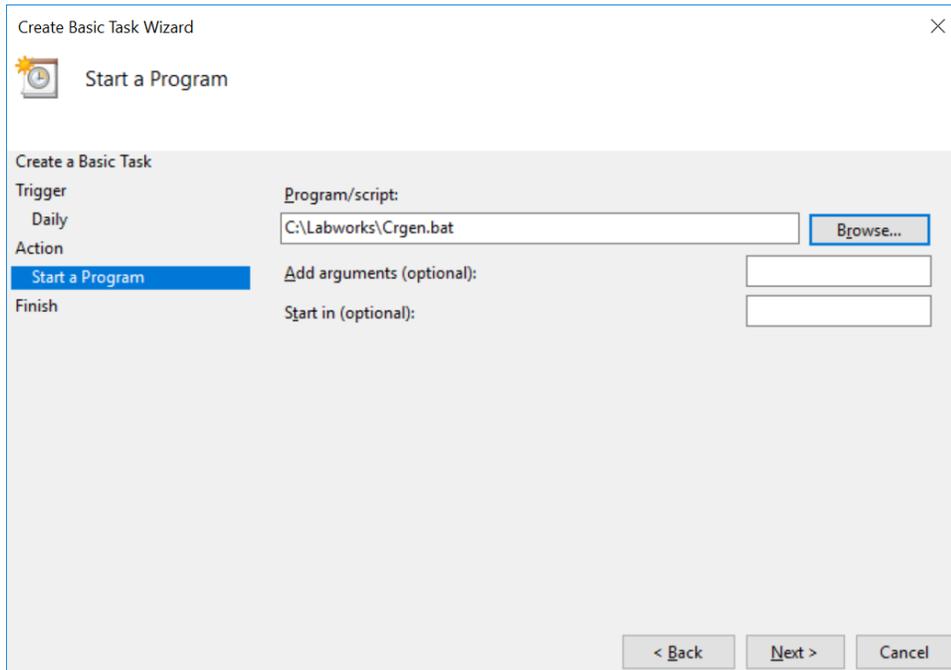

Item Review, Check Open the Properties...

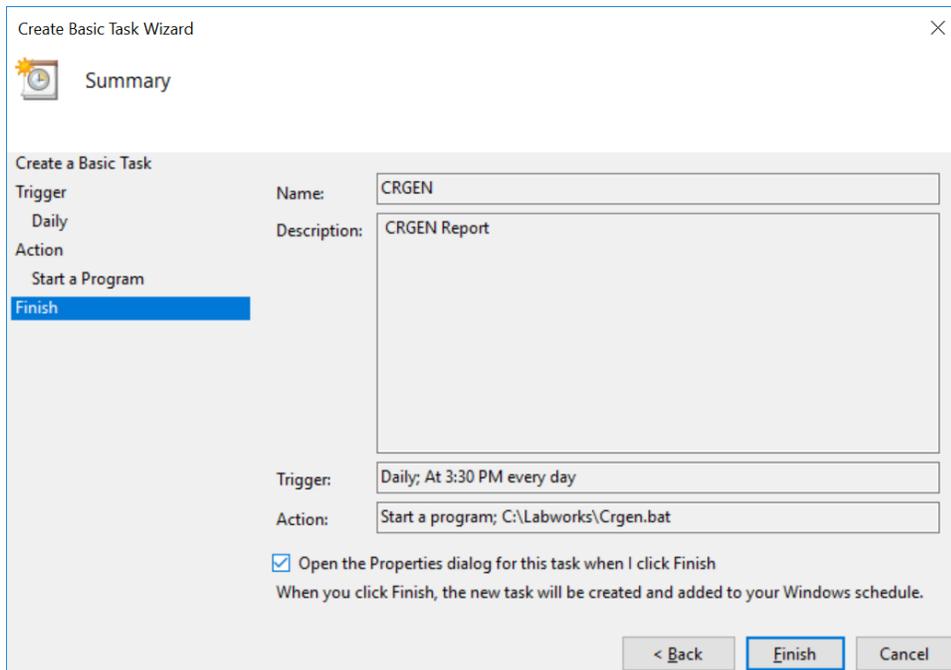

Select "Run whether user is logged on or not"

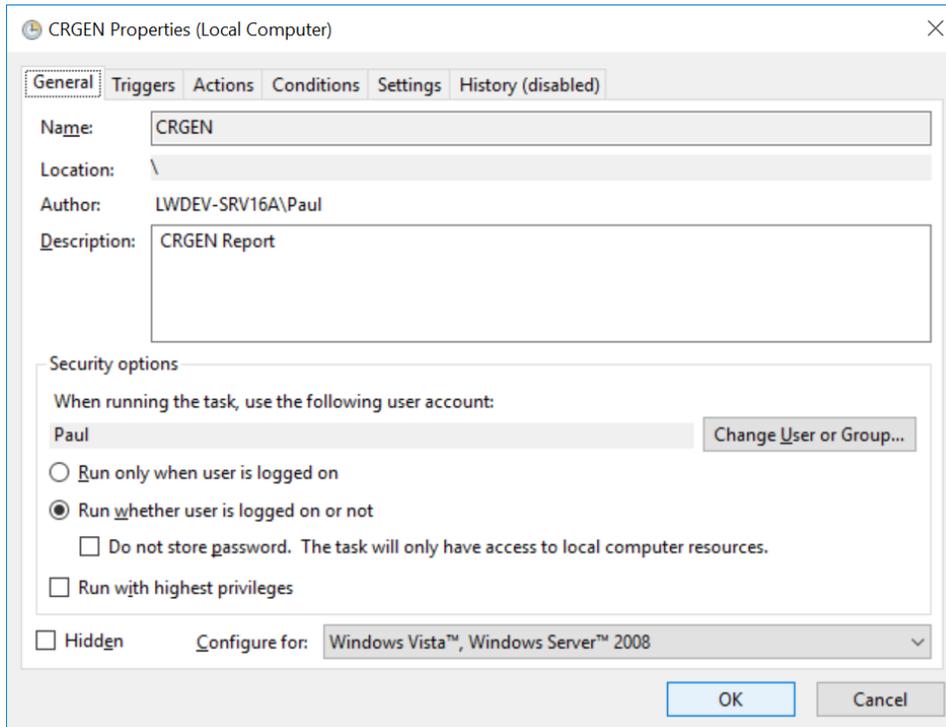

Enter the password for the account when prompted.

