
LABWORKS LIMS v6.8

LABWORKS PICKLISTS

Last Updated: March 30, 2018



LABWORKS PICK-LISTS

1) Default Pick lists	3
2) Reserve Pick list names	4
3) Manage Pick list.....	5

1) Default Pick lists

Pick list which are added into the Database using XSP file.

1) Pick list name: REPORTTYPE

PLITEMNO	PLITEM	DISCRIPTION
0	REPORTTYPE	Report type picklist
1	RECEIPTS	RECEIPTS
2	ROUTESHEETS	ROUTESHEETS
3	WORKSHEETS	WORKSHEETS
4	REPORTTEMPLATE	REPORTTEMPLATE

Used in Applications: LABWORKS Reporting

2) Pick list name: BAUDRATELIST

PLITEMNO	PLITEM	DISCRIPTION
0	Baud Rate List	
1	300	
2	600	
3	1200	
4	2400	
5	4800	
6	9600	

Used in Applications: LABWORKS Instrument configuration

3) Pick list name: DATABITLIST

PLITEMNO	PLITEM	DISCRIPTION
0	Data Bit List	
1	7	
2	8	

Used in Applications: LABWORKS Instrument configuration

4) Pick list name: STOPBITLIST

PLITEMNO	PLITEM	DISCRIPTION
0	Stop Bit List	
1	One	
2	OnePointFive	
3	Two	

Used in Applications: LABWORKS Instrument configuration

5) **Pick list name:** RLTCODES

PLITEMNO	PLITEM	DISCRIPTION
0	RLT Codes	
1	ND	Not detected

Used in Applications: LWRoundOffEngine, LWWebTop

2) [Reserve Pick list names](#)

Following are the optional picklists which are used in different LABWORKS applications.

1) **Pick list name:** PICKLIST_EXC

Used in Applications: LABWORKS Analysis Maintenance

2) **Pick list name:** PICKLIST_INC

Used in Applications: LABWORKS Analysis Maintenance

3) **Pick list name:** SPECIALIST NO.

Used in Applications: FLAG WindowsSMS (Custom application)

4) **Pick list name:** PSDEPT

Used in Applications: LWSampleLogin, Process Scheduler

5) **Pick list name:** SC_QA_PREFIX **Used in**

Applications: NGIIAdapter

6) **Pick list name:** C_QA_SUFFIX **Used in**

Applications: NGIIAdapter

7) **Pick list name:** ITEM_UNITS

Used in Applications: LWChemicalInventory

8) **Pick list name:** SHEFLIFE_UNIT **Used in**

Applications: LWChemicalInventory

9) **Pick list name:** ITEM_TYPE

Used in Applications: LWChemicalInventory

10) **Pick list name:** SLIM_LABS

Used in Applications: LABWORKS SLIMLINK (Custom application)

11) **Pick list name:** MRB_COURIER_STATUS

Used in Applications: MRBE_BILLING (Custom application)

12) **Pick list name:** SAMPNOTCOLREASONS

Used in Applications: LABWORKS Sample Collection App


13) **Pick list name:** USER_SITES

Used in Applications: LWLicenseManagement


3) Manage Pick list

a) Create new Pick list -

- Open Picklist Maintenance application.

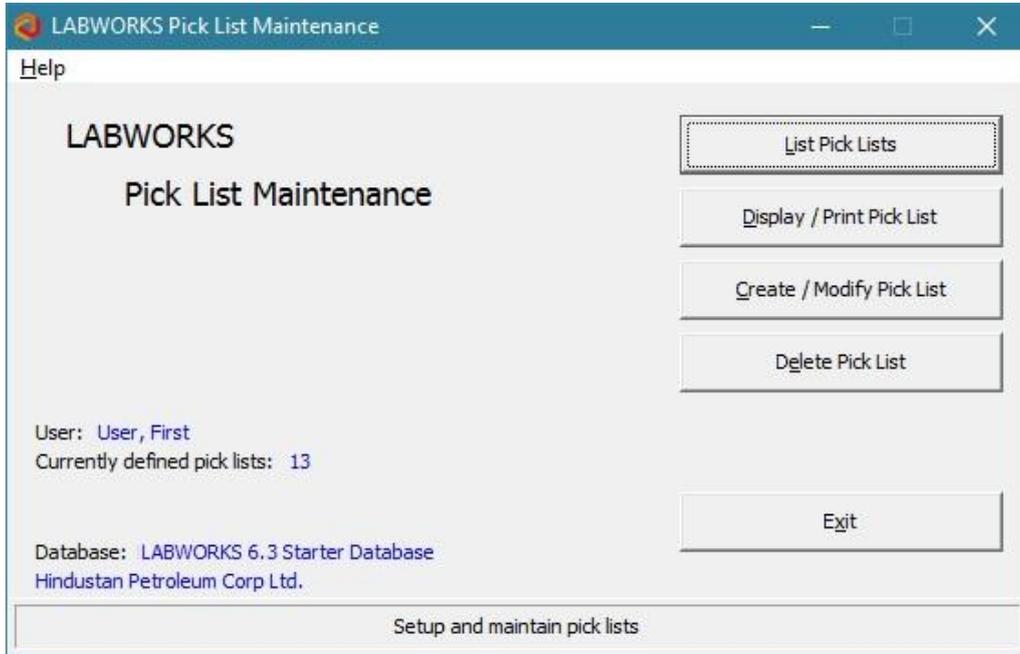


- Click on Create/Modify Pick List. Below window will appear.

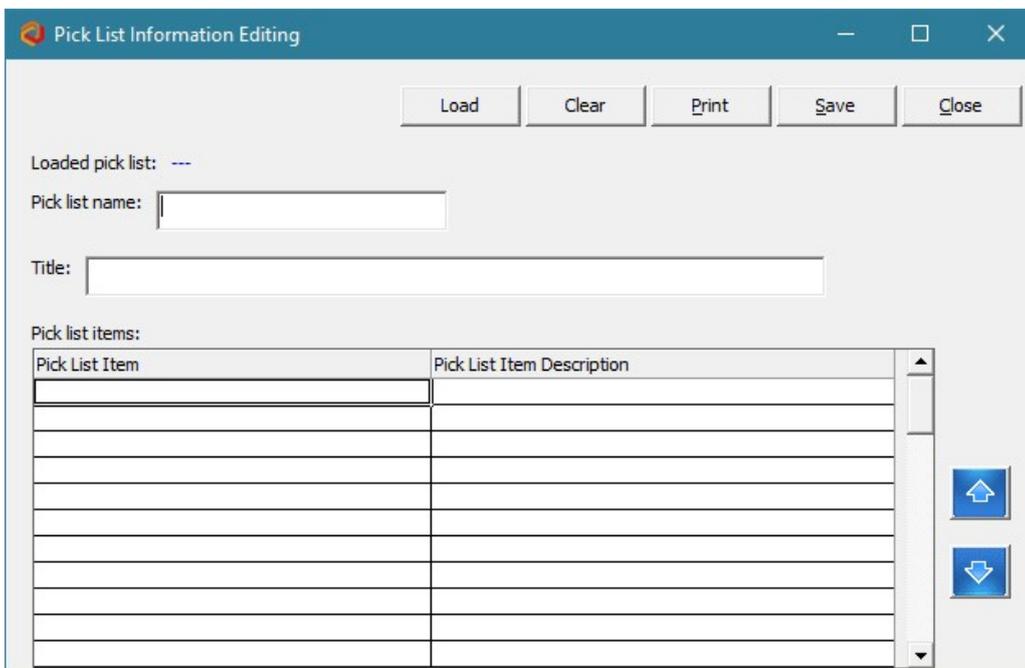


Enter Pick list name, title and picklist data in the “Pick list items” grid and click on “Save” button to save a pick list.

b) Edit Pick list –


- Open Picklist Maintenance application.

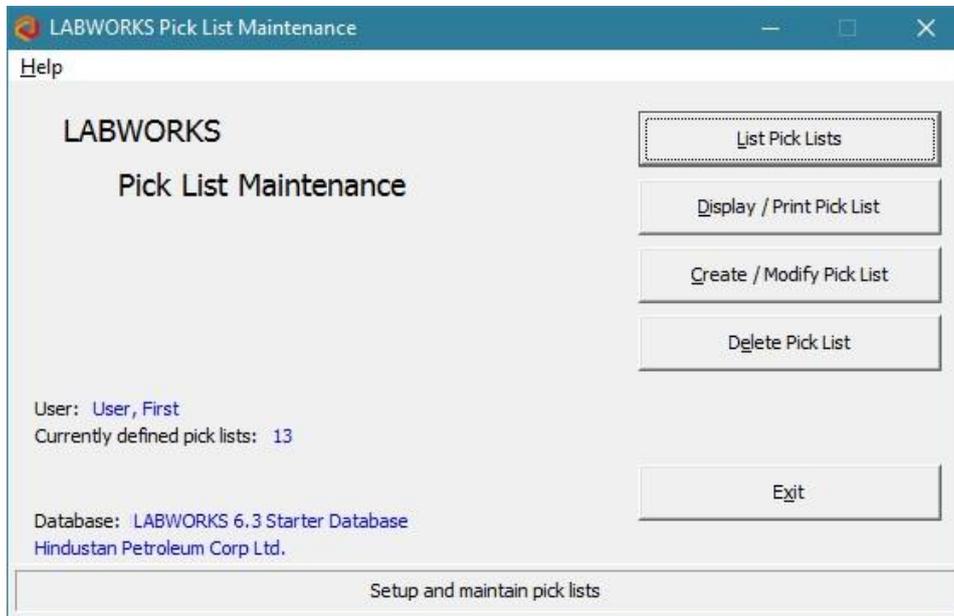


- Click on Create/Modify Pick List. Below window will appear.

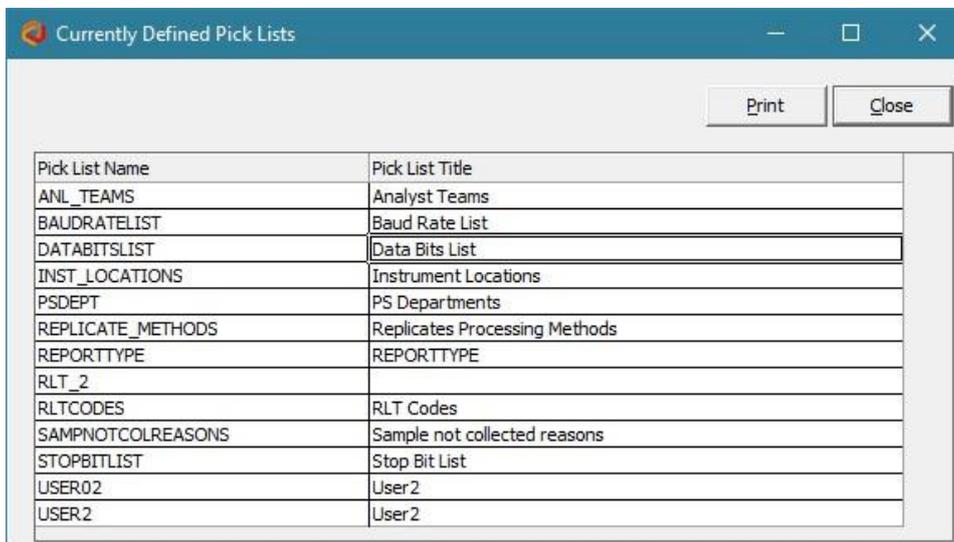

- Click on “Load” button, below window will appear.

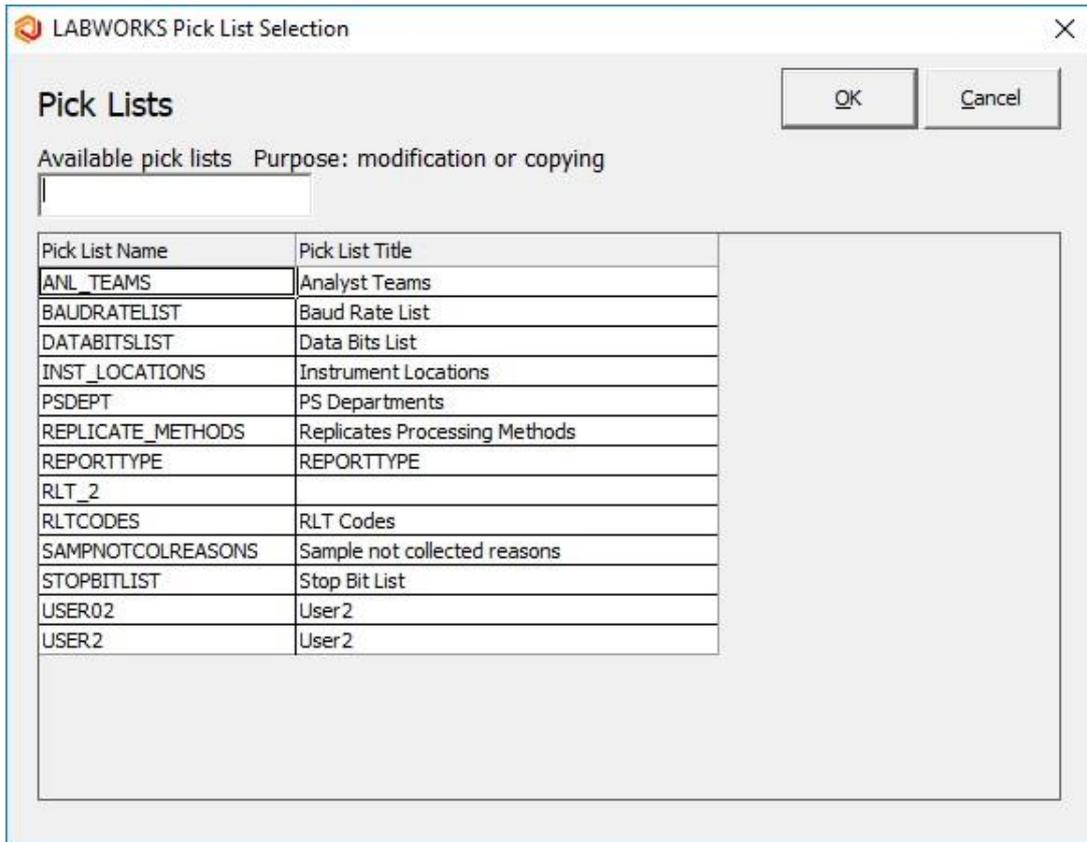


- Select picklist to edit and click “OK”. Selected Pick list will be open in edit mode.

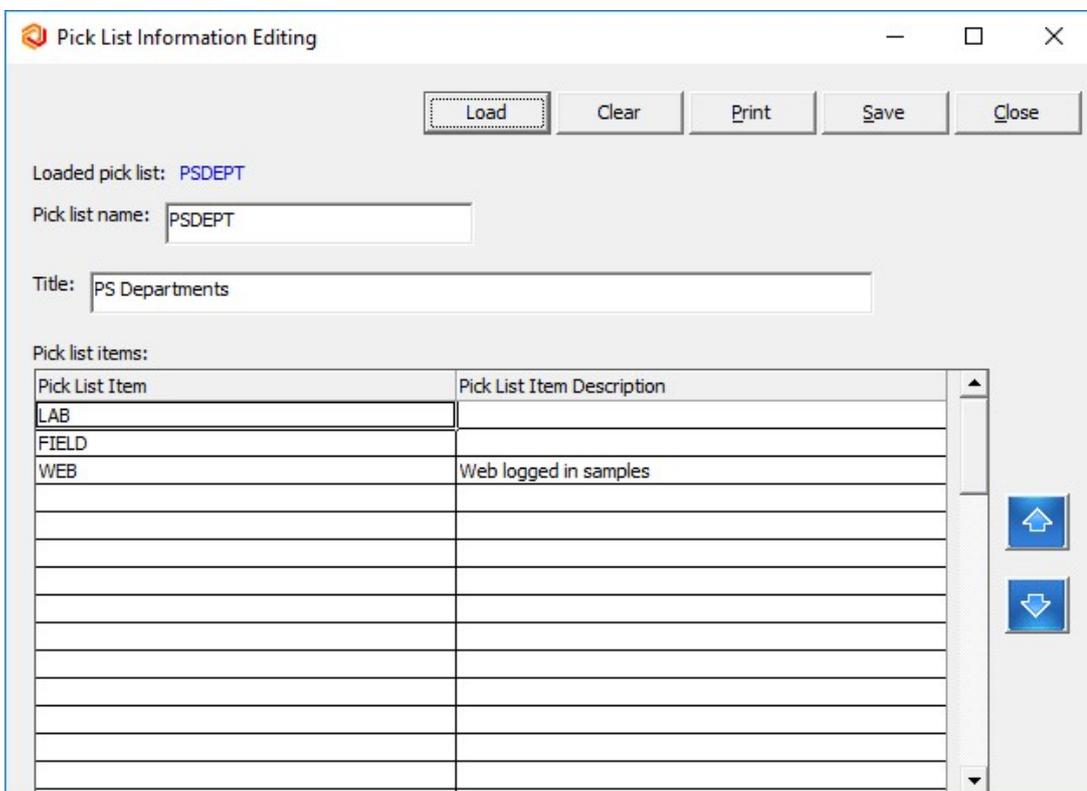


After modification, click on “Save” button to update the Pick list.

c) Delete Pick list –


- Open Picklist Maintenance application.

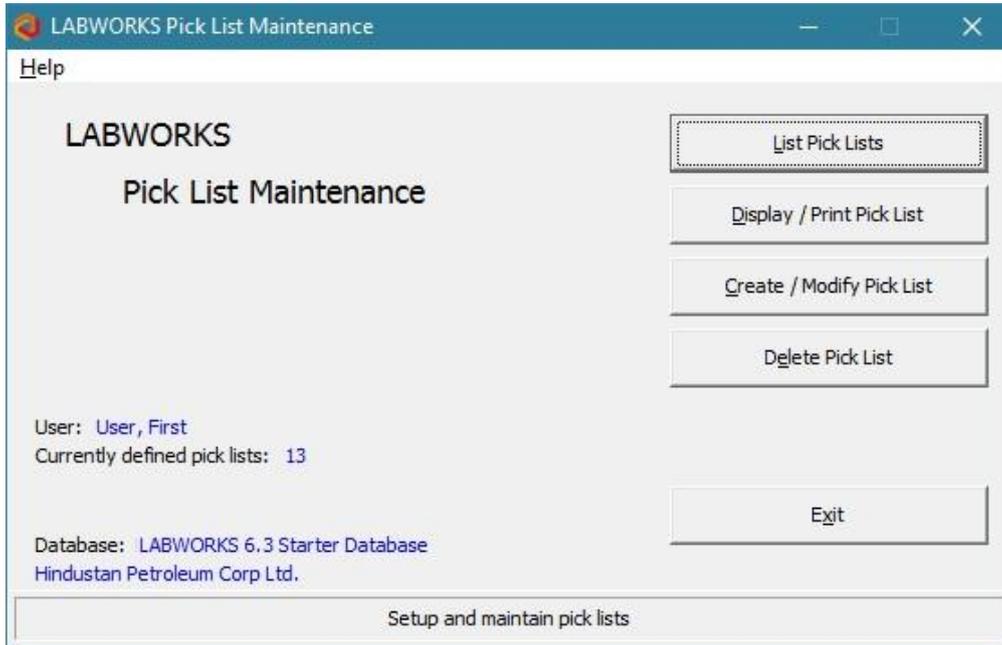

- Click on “Delete Pick List” button, below window will appear.

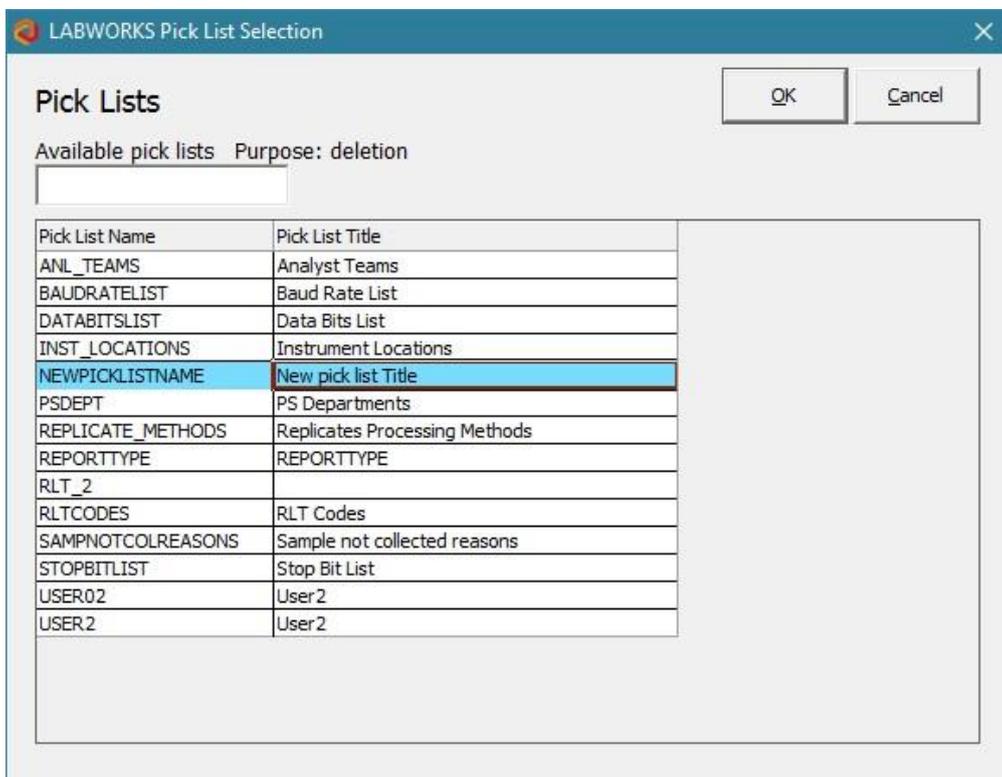

- Select pick list to delete and click OK

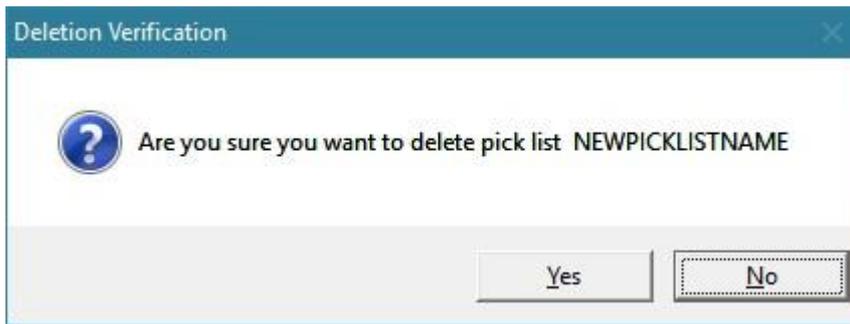

A confirmation window will appear. Click on "Yes" to delete Pick list permanently.